EN

50 51 T probe

Refer to the meter manual for more information.

Caution

⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*

⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

Introduction

This probe is a combination gel-filled electrode with a built-in temperature sensor for pH measurements in wastewater, foods and samples with high solid content.

Technical information

Range	2 to 14 pH
Junction	Open
Reference	Ag/AgCl
Temperature	Pt 1000
Minimum depth	15 mm (0.6 in.)
Operating temp	0–80 °C (32–176 °F)
Limits	Low-conductivity or pH < 2
Connector type	MP-5

Preparation

Install the protective shroud.

Connect the probe

Caution

⚠ *Do not turn or twist the probe connector.*

1 Calibration

Caution

⚠ *Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.*

⚠ *Rinse sensor thoroughly between standards.*

For 1 or 2-point pH calibration: push ✓ after 1 or 2 standards.

2 Sample measurement

Continuous: push ✓ 2 times.

Change the parameter: push and hold ✓ during stabilization.

Maintenance and storage

Cleaning: soak the probe in a cleaning solution. For proteins, soak in 2% NaOH for 20 min. For fats and oils, wash with ethanol or a soap solution at 40 °C. Rinse with deionized water.

Storage: put the probe in the protector with KCl 3 M solution.

Troubleshooting

Be sure that the correct probe type is used. Do not hold the tubes by hand to prevent temperature changes.

Warranty

6 months for manufacturing faults only. Damage from use is not covered.

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