50 55

User Manual

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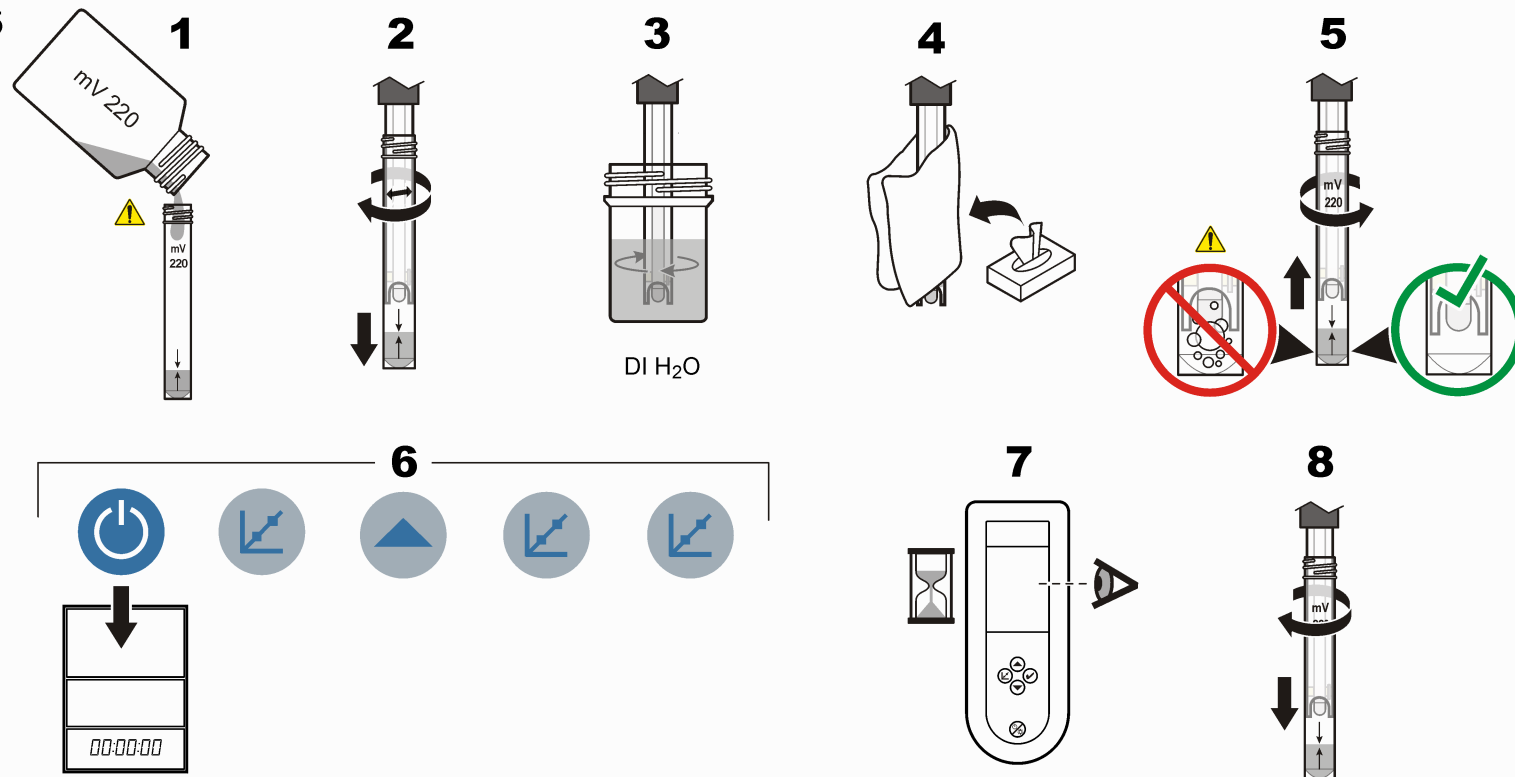
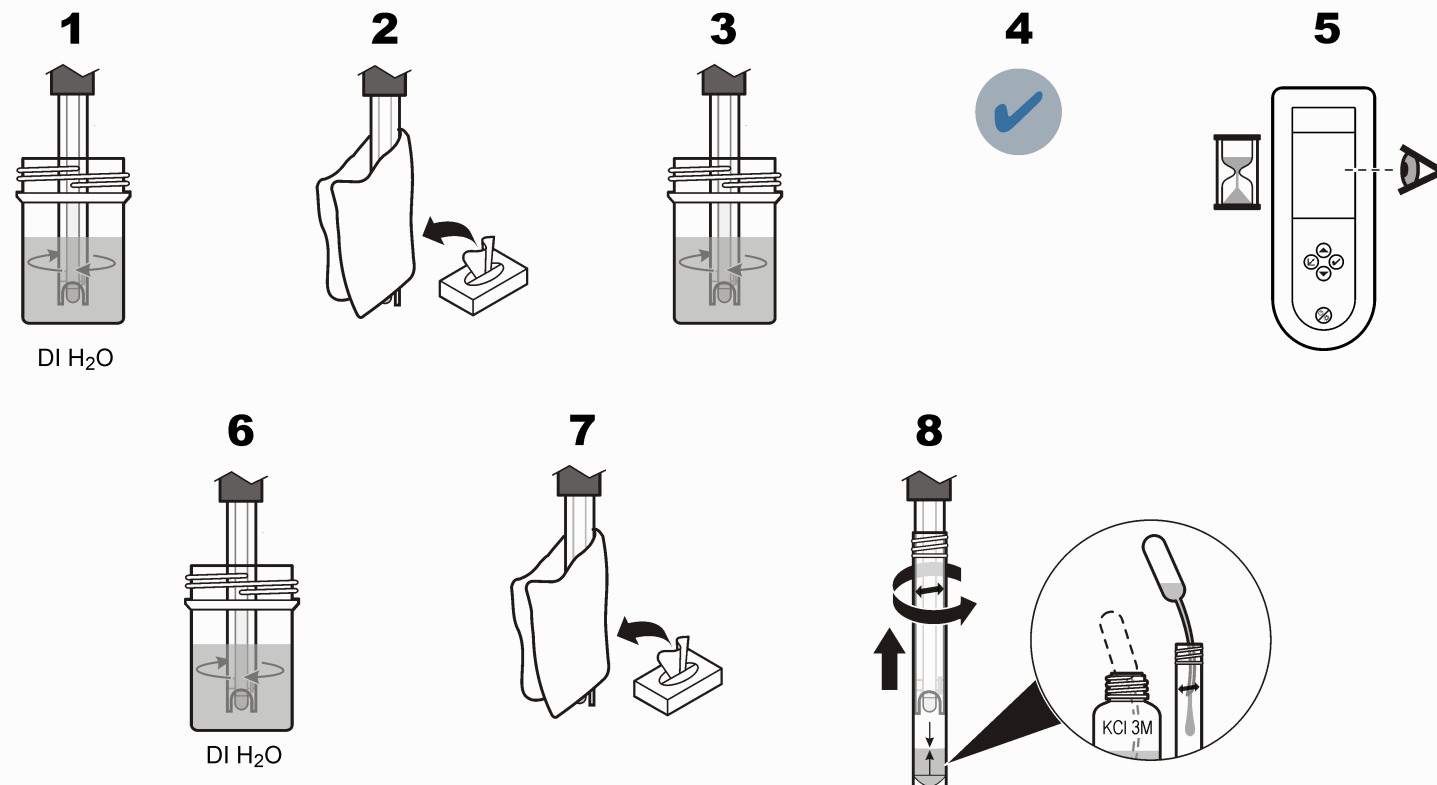
50 55

**2****50 55 probe**

Refer to the meter manual for more information.

Caution

⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*

⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

Introduction

This probe is a combination gel-filled electrode for ORP measurements in general aqueous samples, wastewater treatment and galvanic baths.

Technical information

Range	± 2000 mV
Junction	Ceramic
Reference	Ag/AgCl
ORP electrode	Platinum
Minimum depth	18 mm (0.7 in.)
Operating temp	0 to 80 °C (32 to 176 °F)
Limits	Low-conductivity, dirty or viscous solutions
Connector type	MP-5

Connect the probe**Caution**

⚠ *Do not turn or twist the probe connector.*

1 Calibration**Caution**

⚠ *Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.*

2 Measurement

Continuous: push ✓ 2 times.

Change the parameter: push and hold ✓ during stabilization.

Maintenance and storage

Cleaning: soak the probe in a cleaning solution. Rub the surface of the platinum electrode with a soft abrasive, for example toothpaste. Rinse with deionized water.

Storage: put the probe in the protector with KCl 3 M solution.

Troubleshooting

Be sure that the correct probe type is used. Do not hold the tubes by hand to prevent temperature changes.

Warranty

6 months for manufacturing faults only. Damage from use is not covered.

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