



Intro to **EnSURE™** Touch

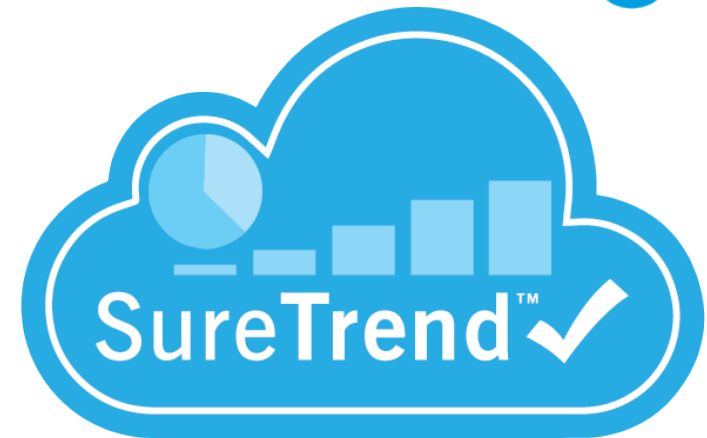

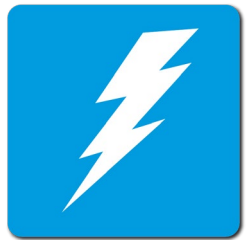


Quick Test

- Run a test in as few as two touches








Quick Test

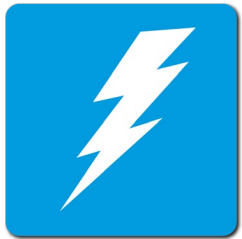

Quick Test

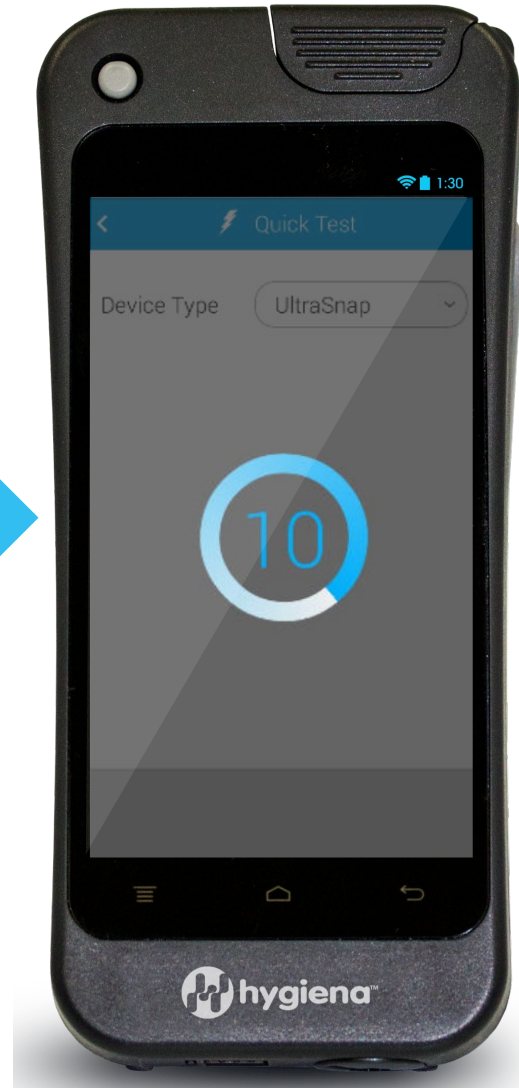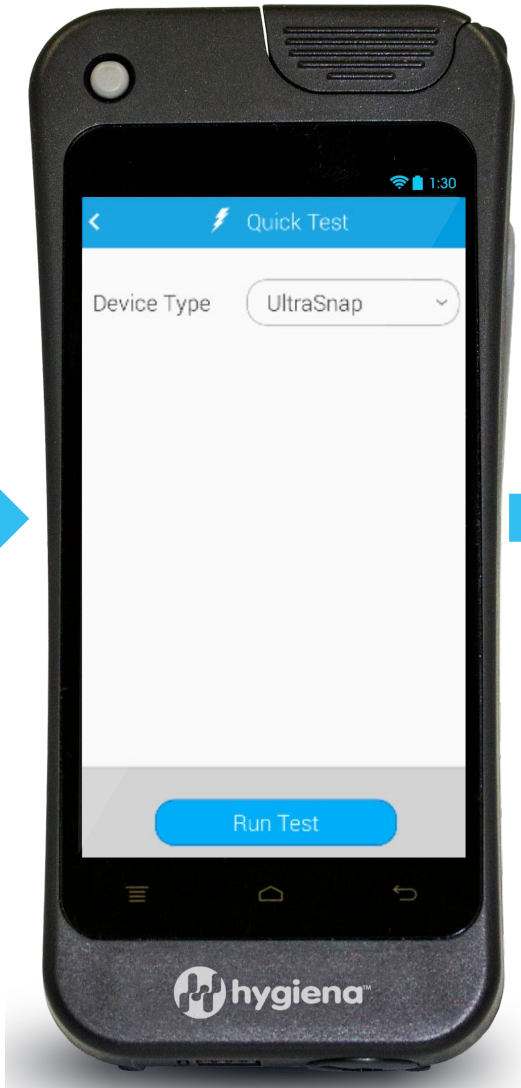

EnSURE™ Touch

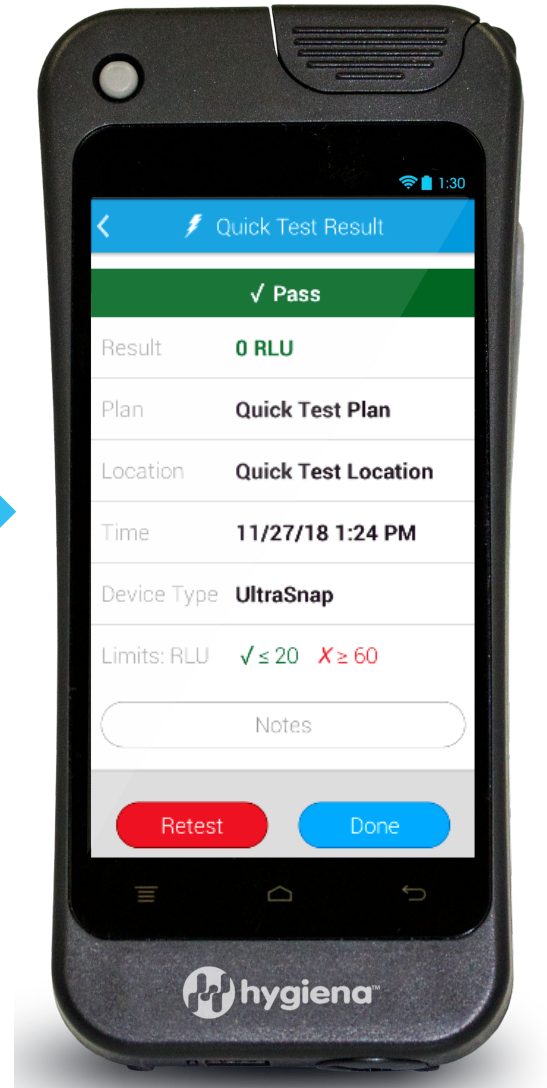



Users

- Manage user capabilities with roles
 - Operators perform testing
 - Administrators have full access and can perform testing
- Password protection








Adding Users

EnSURE™ Touch



Users

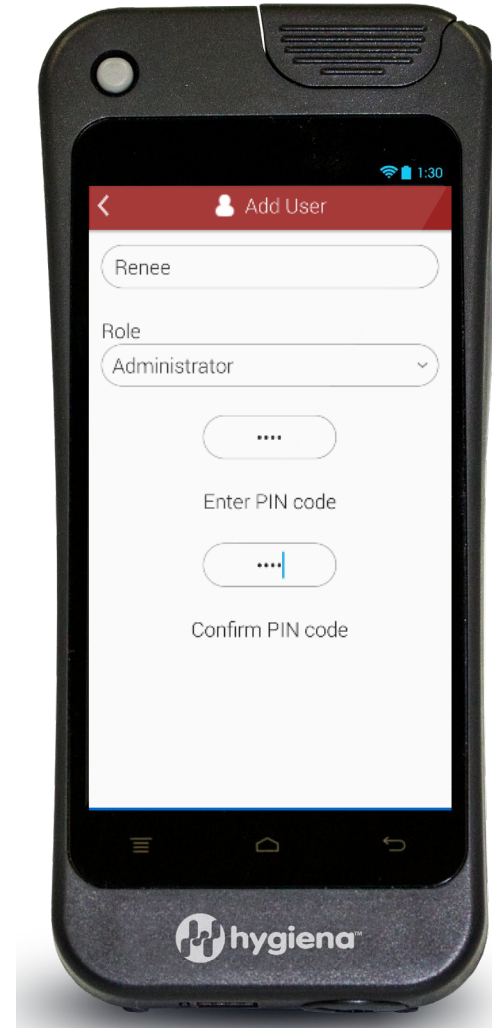
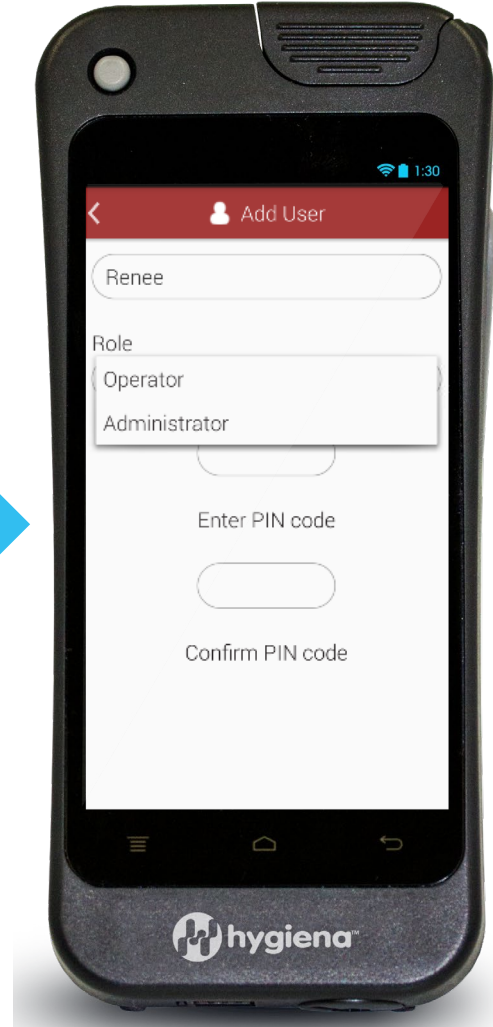
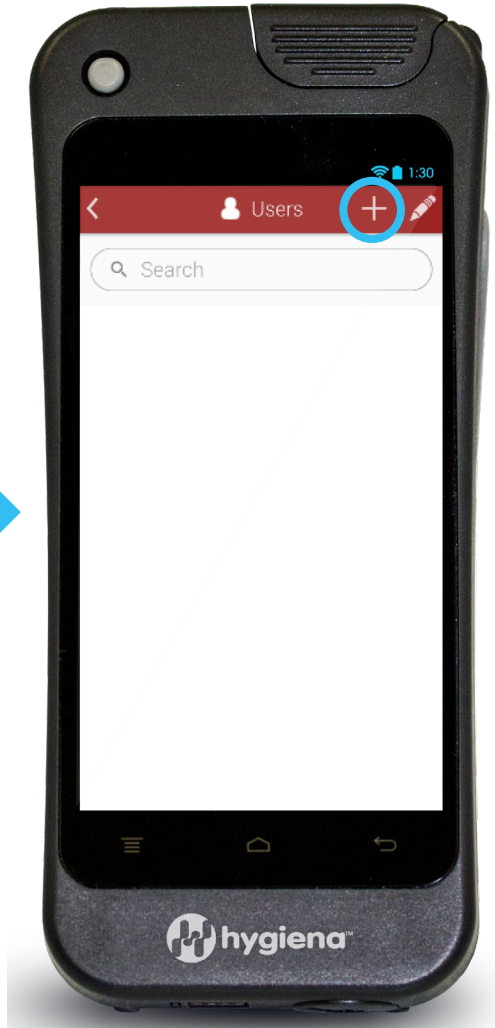




Managing Users



Users

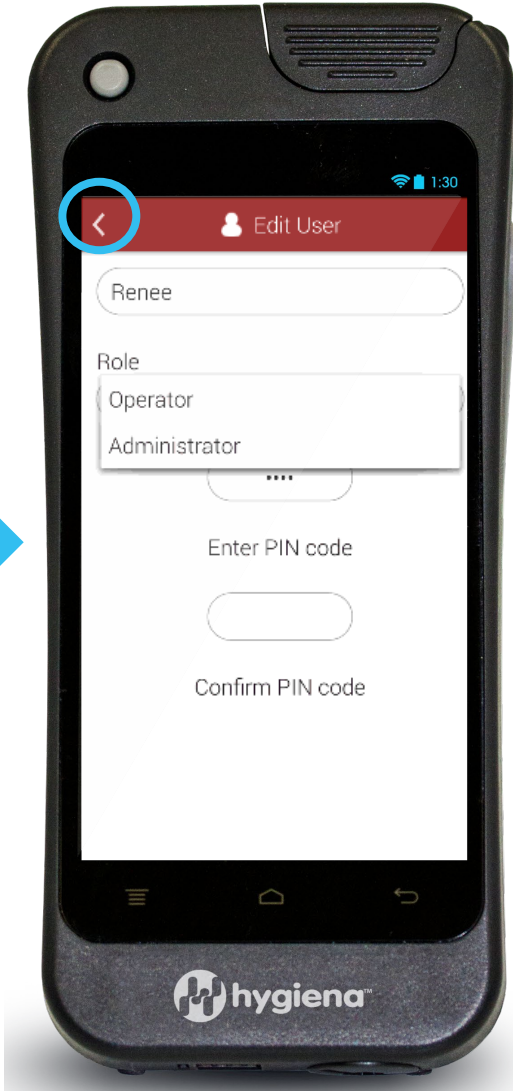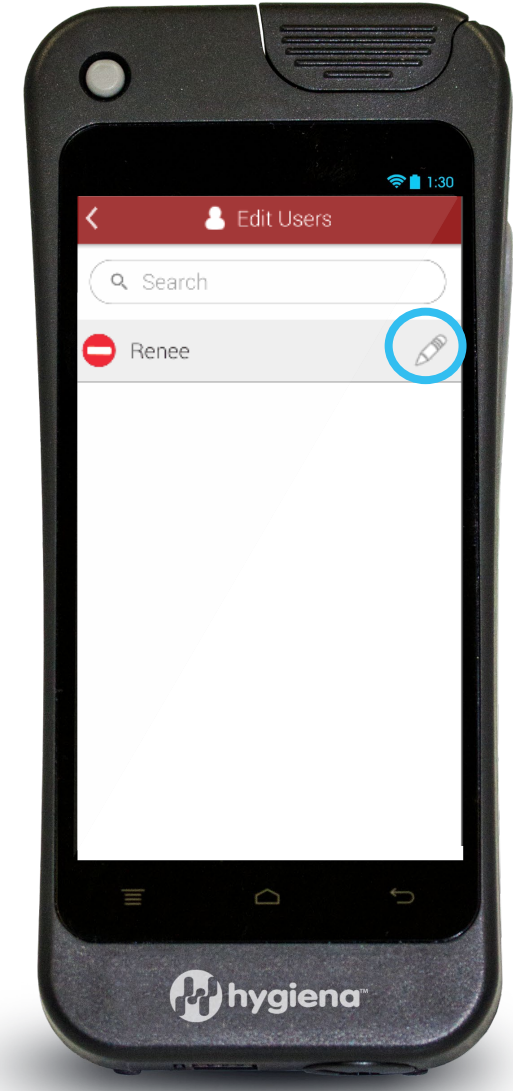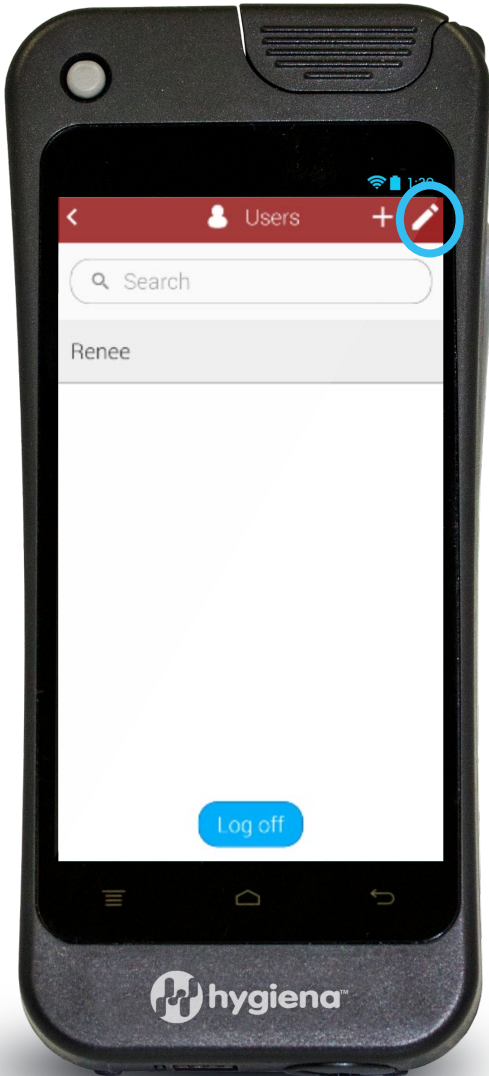

Don't worry, pressing the back button saves your work

EnSURE™ Touch



Plans

- Create plans directly on the instrument without needing to go to your computer to use software
- Randomize testing locations in the plan to ensure sample coverage
- Schedule recurring plans








Creating a Plan



Plans


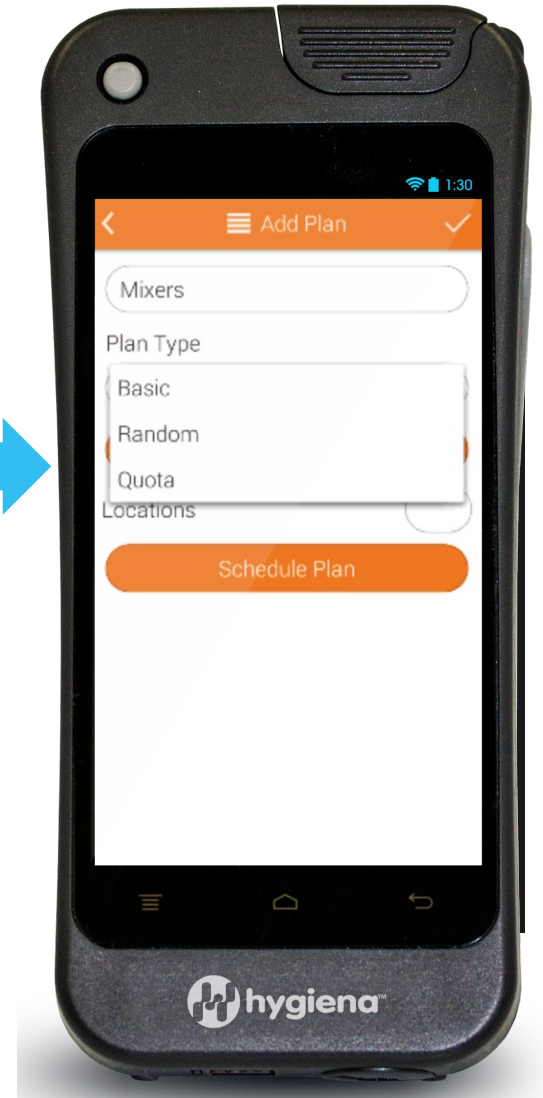
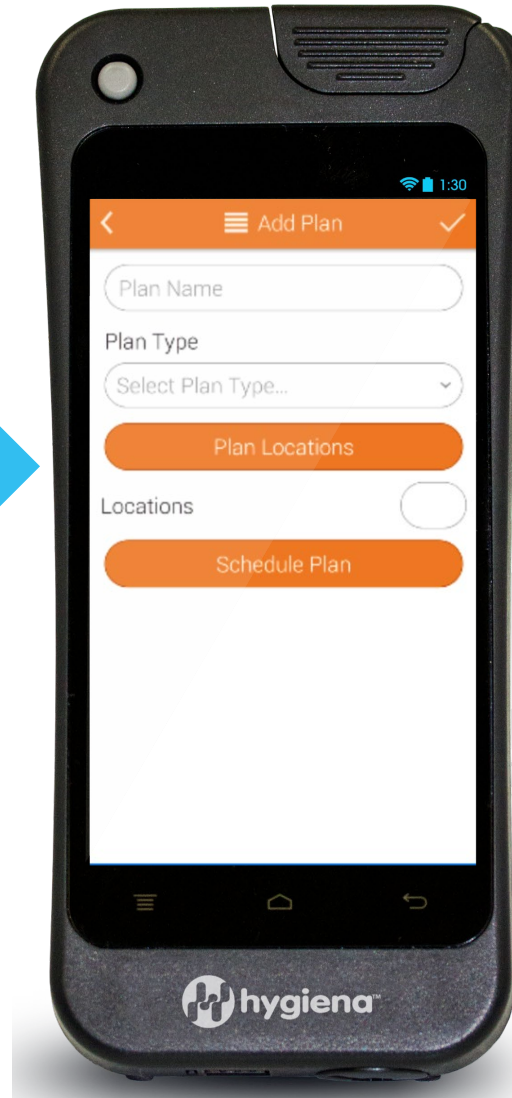
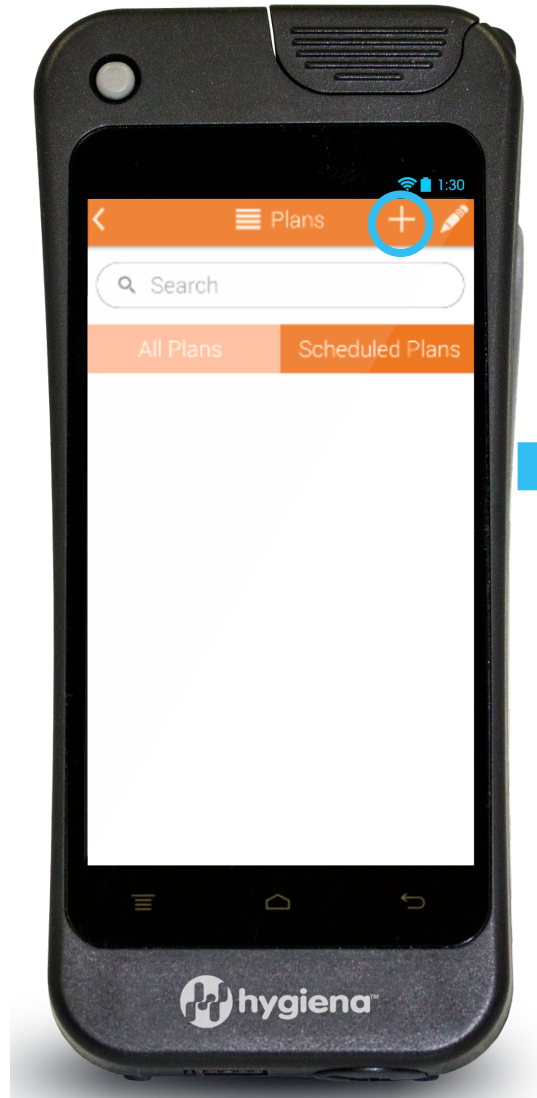


EnSURE™ Touch



Adding Locations to a Plan

EnSURE™ Touch

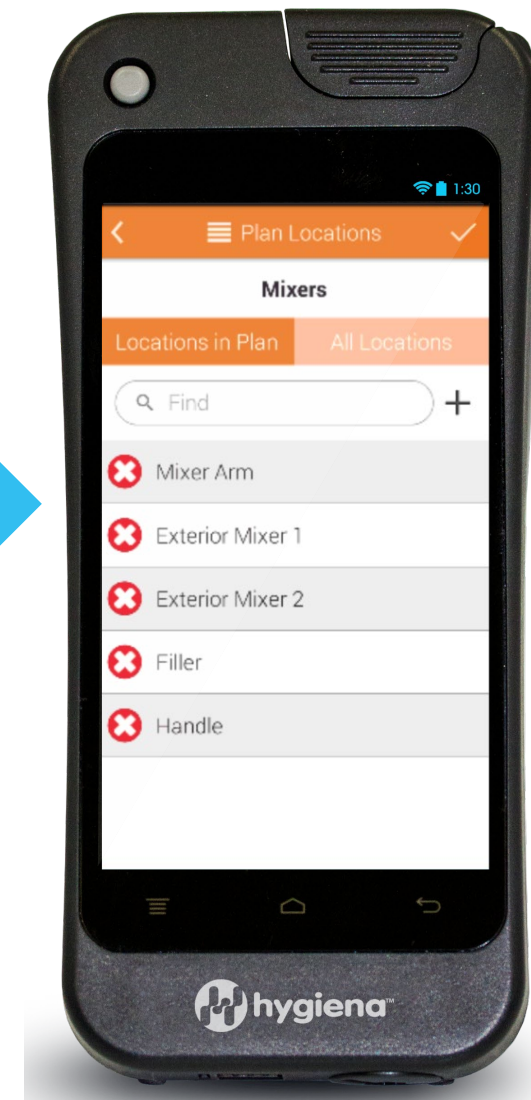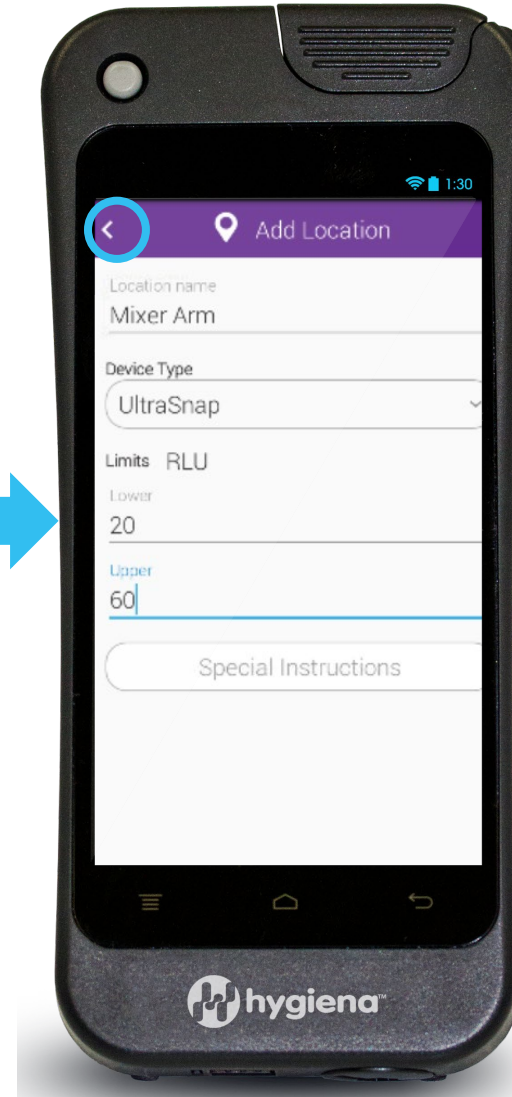
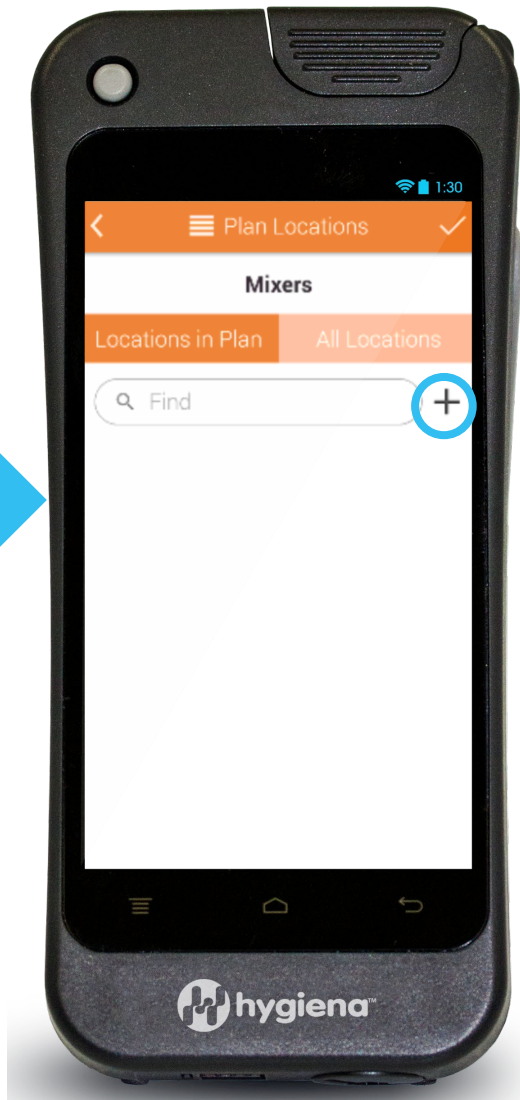
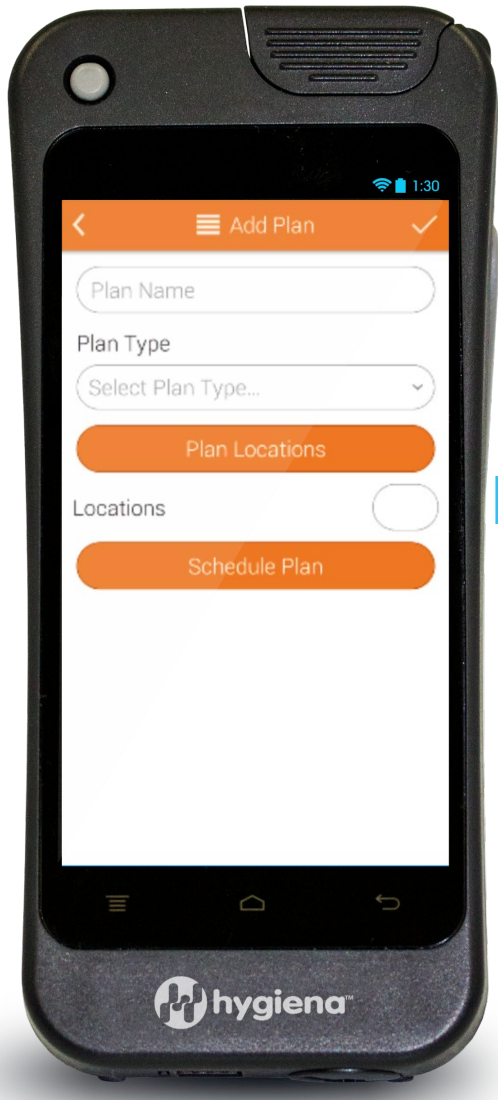


Don't worry, pressing the back button
saves your work



Scheduling a Plan

EnSURE™ Touch

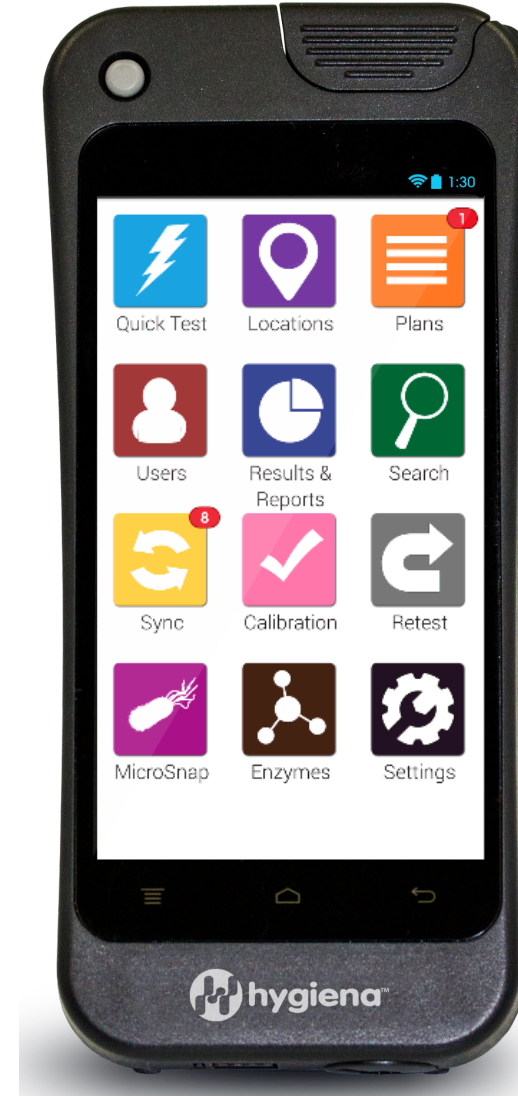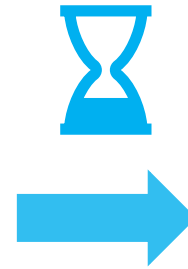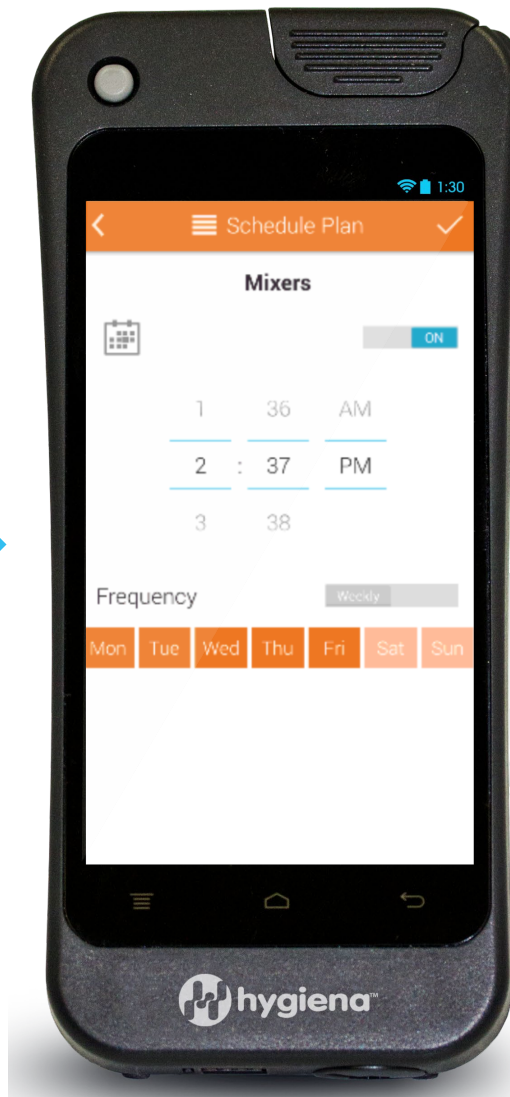

Plans

When it's time to run a scheduled plan a red icon will appear above the orange plan icon






Retest

- Swipe through & view failed test locations
- Instantly run re-tests
- Easily show effectiveness of corrective action







EnSURE™ Touch

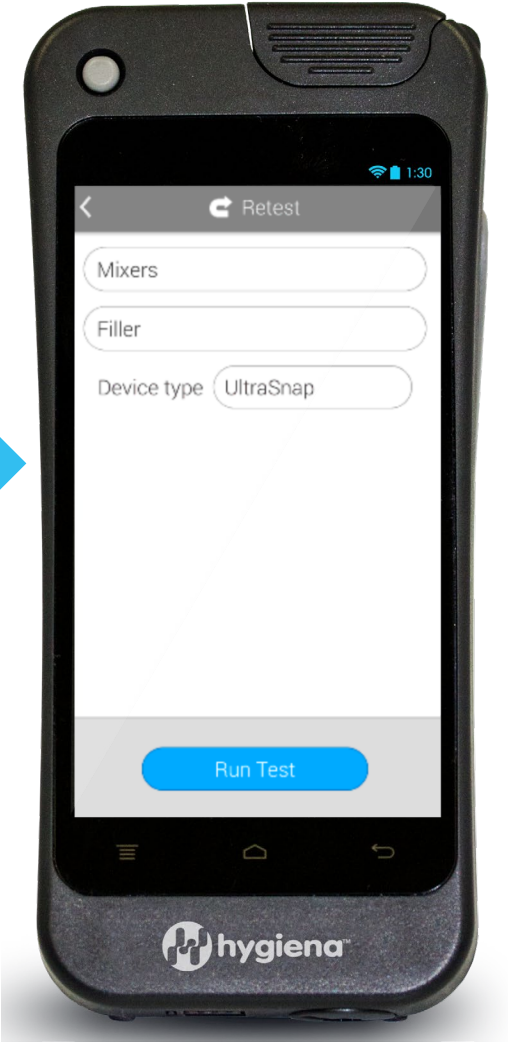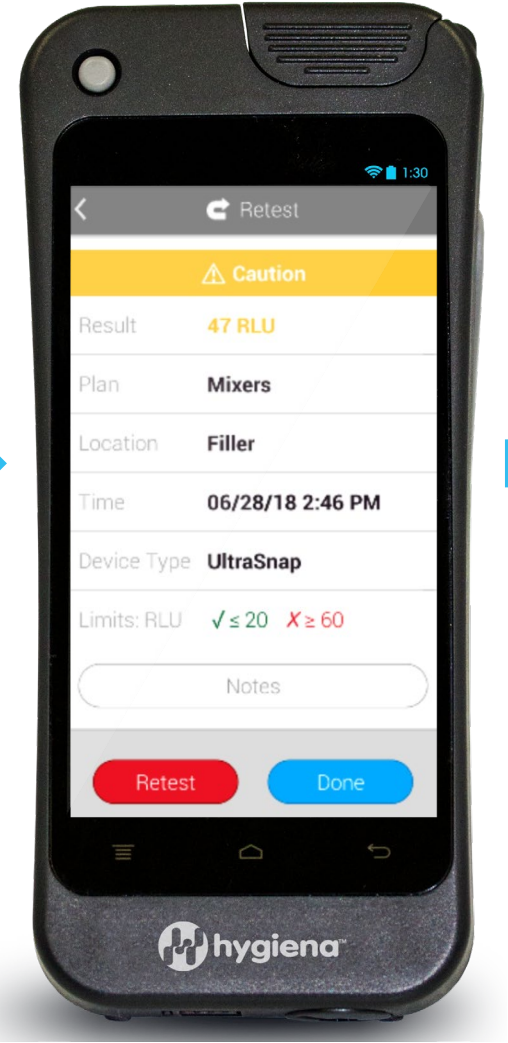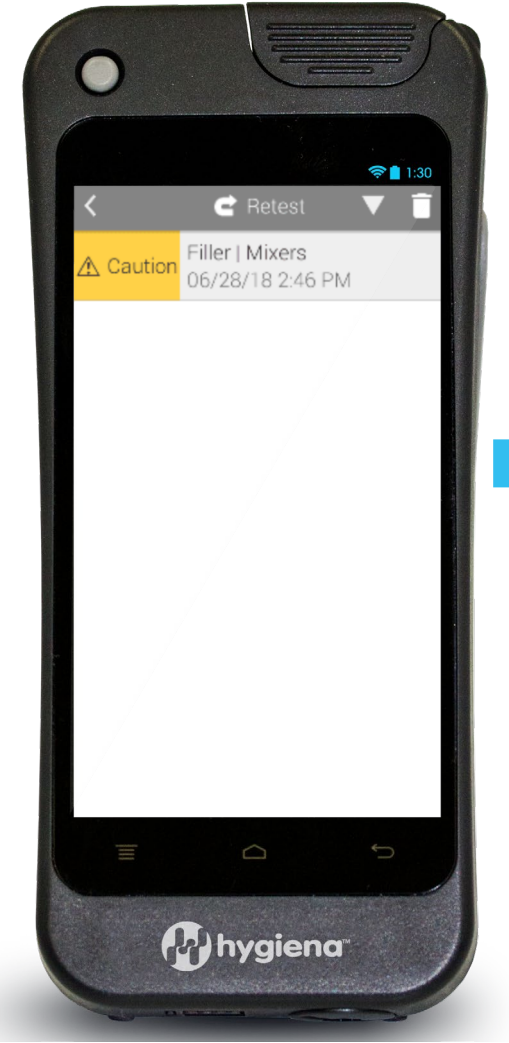



Retest

EnSURE™ Touch

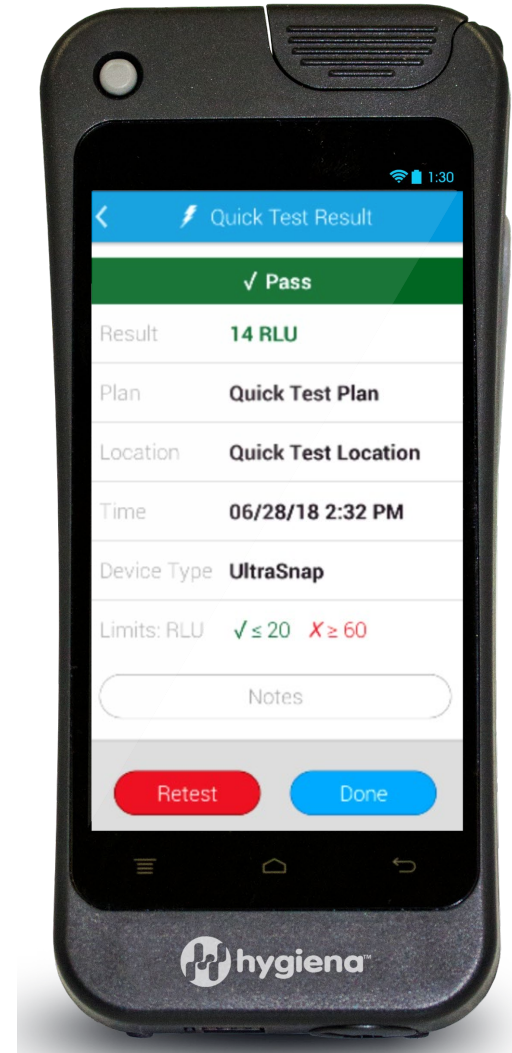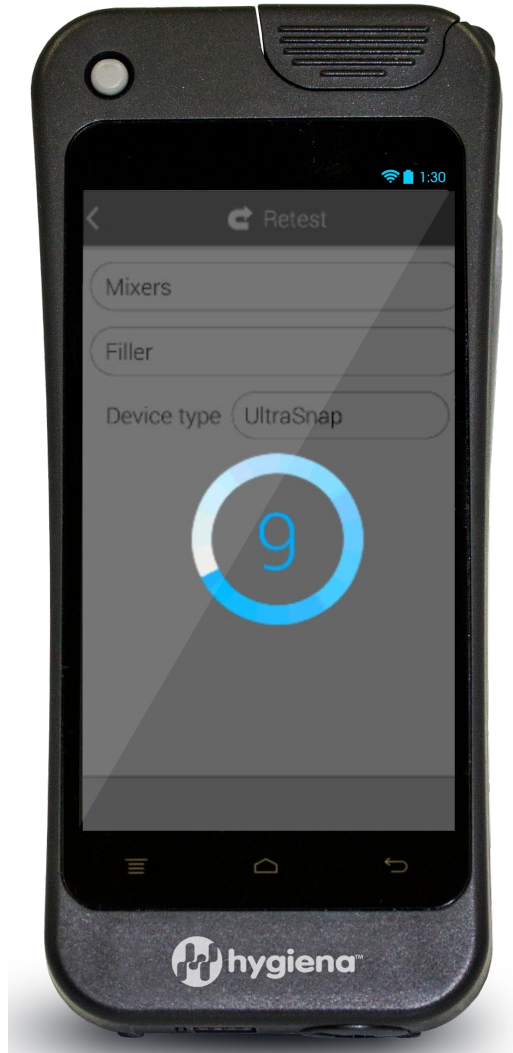



Sync

- Wirelessly sync data to between your EnSURE Touch and SureTrend Cloud account








Turn on Syncing From the EnSURE Touch

EnSURE™ Touch



Settings

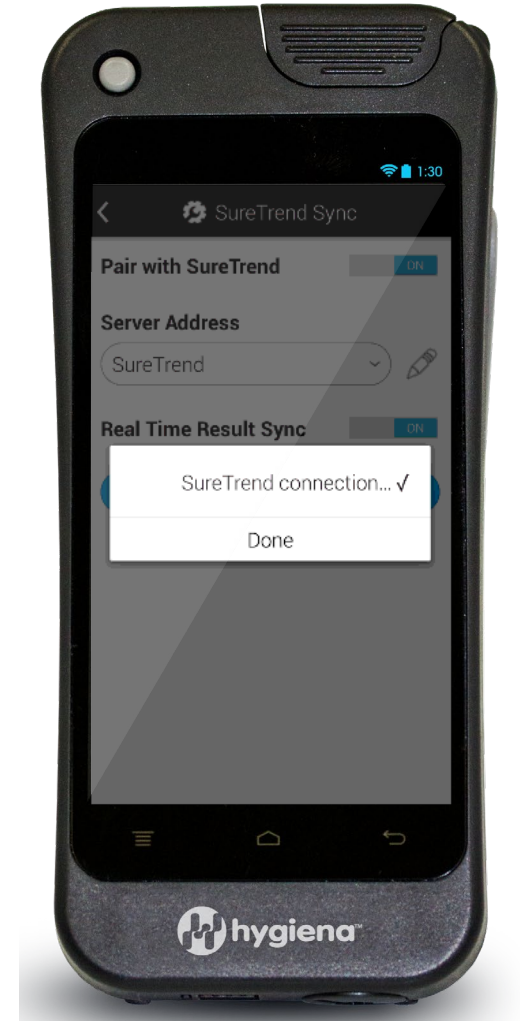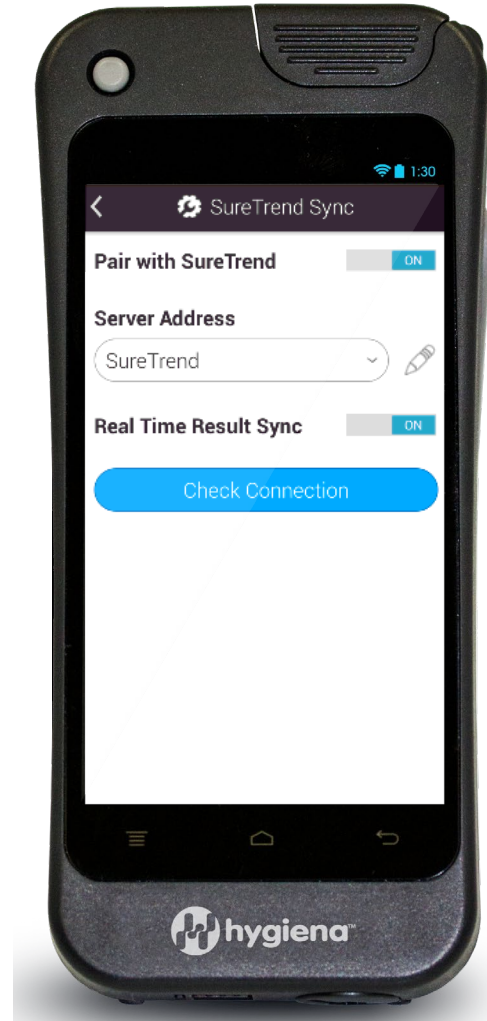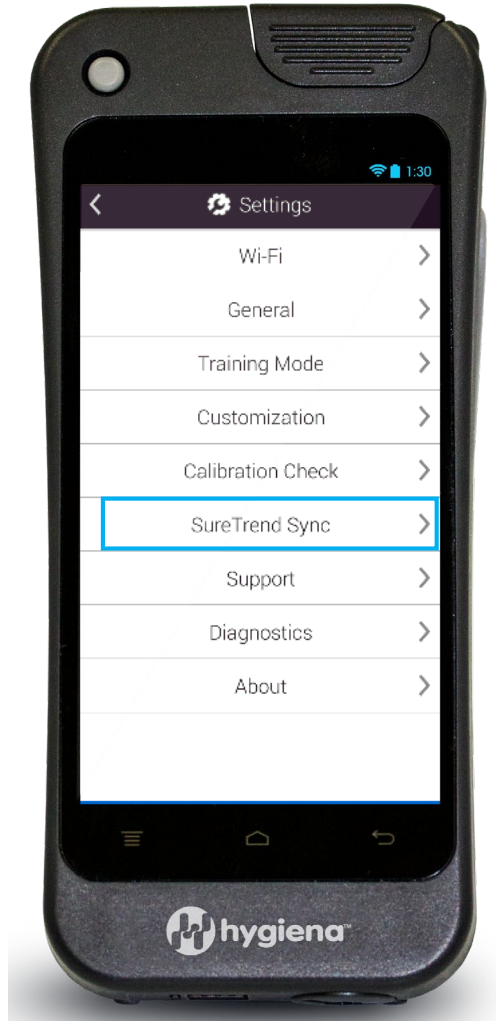

***Instrument must be registered on SureTrend Cloud. Visit help.hygiena.com for tutorial.*






Syncing

EnSURE™ Touch



Sync

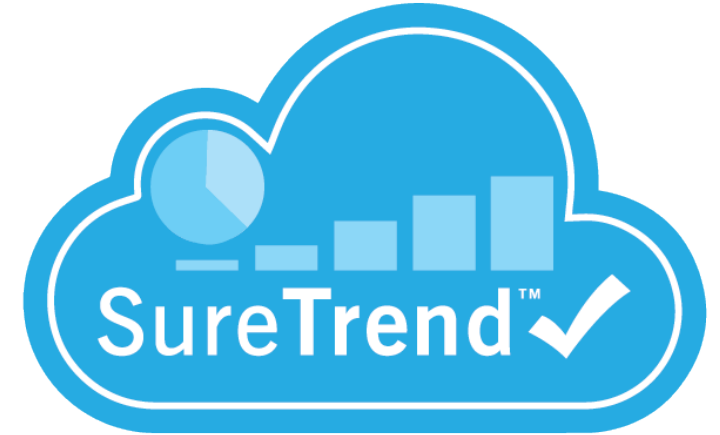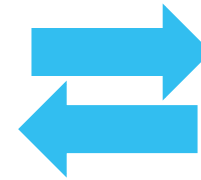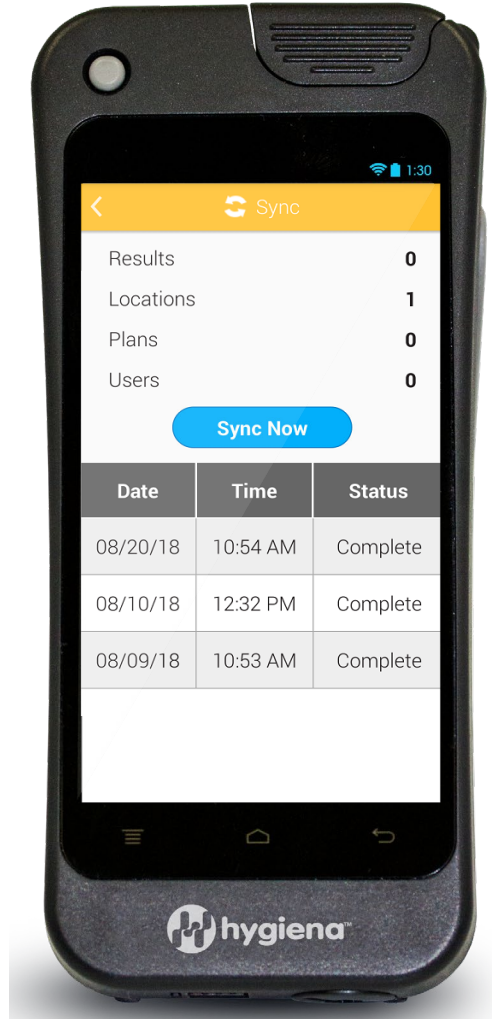

***All results will be synced with SureTrend Cloud. The process should take less than a minute, but may be longer depending on how many results need to be synced.*



Results & Reports

- Basic reporting and data trending available on the instrument for on the go analysis








Results & Reports



Results & Reports

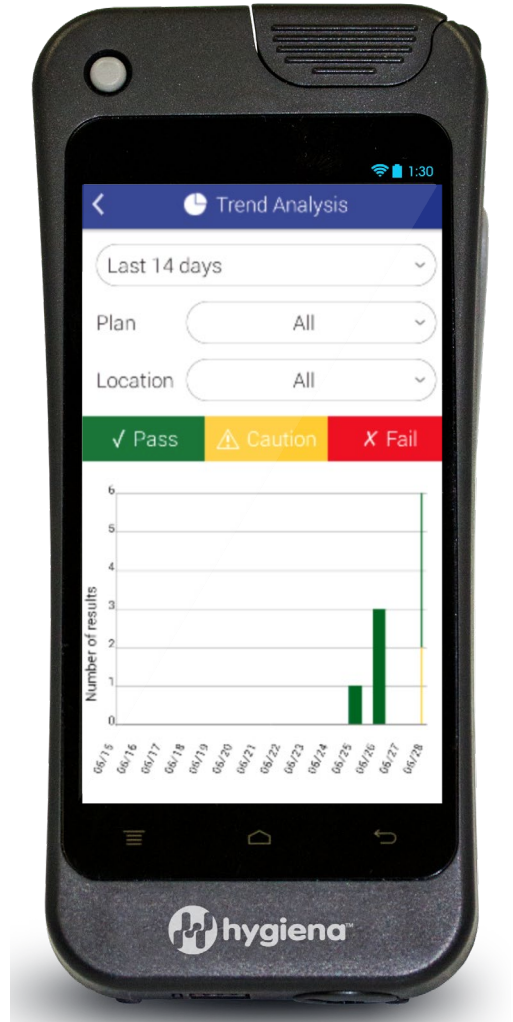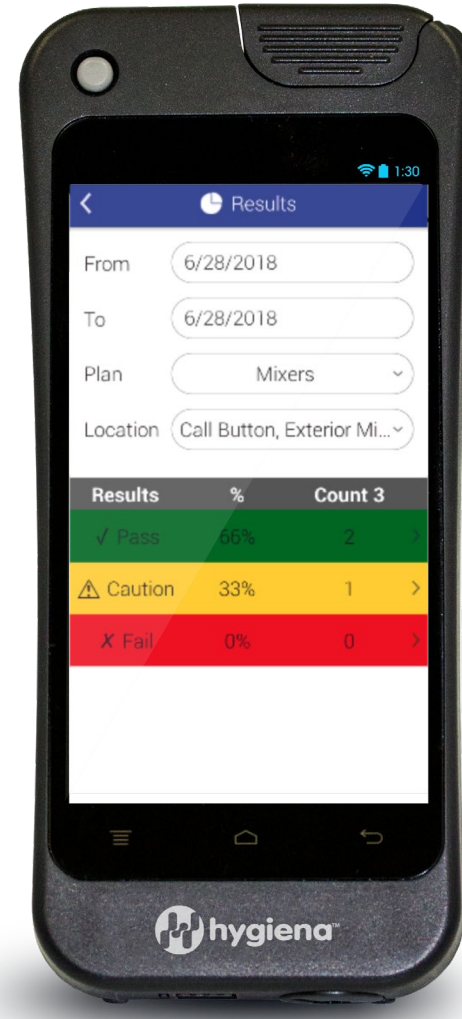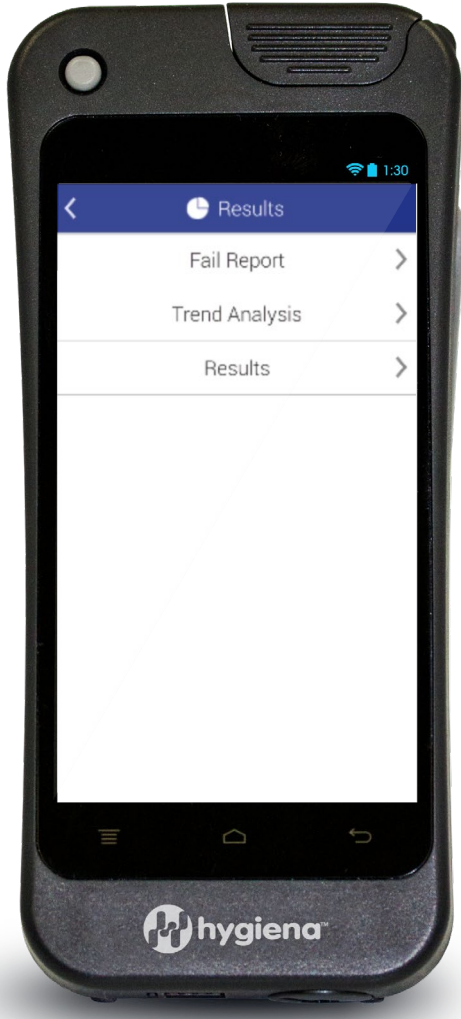

EnSURE™ Touch



RLU Conversion


- Makes it easy to switch to Hygiena without changing upper and lower threshold limits







RLU Conversion

EnSURE™ Touch

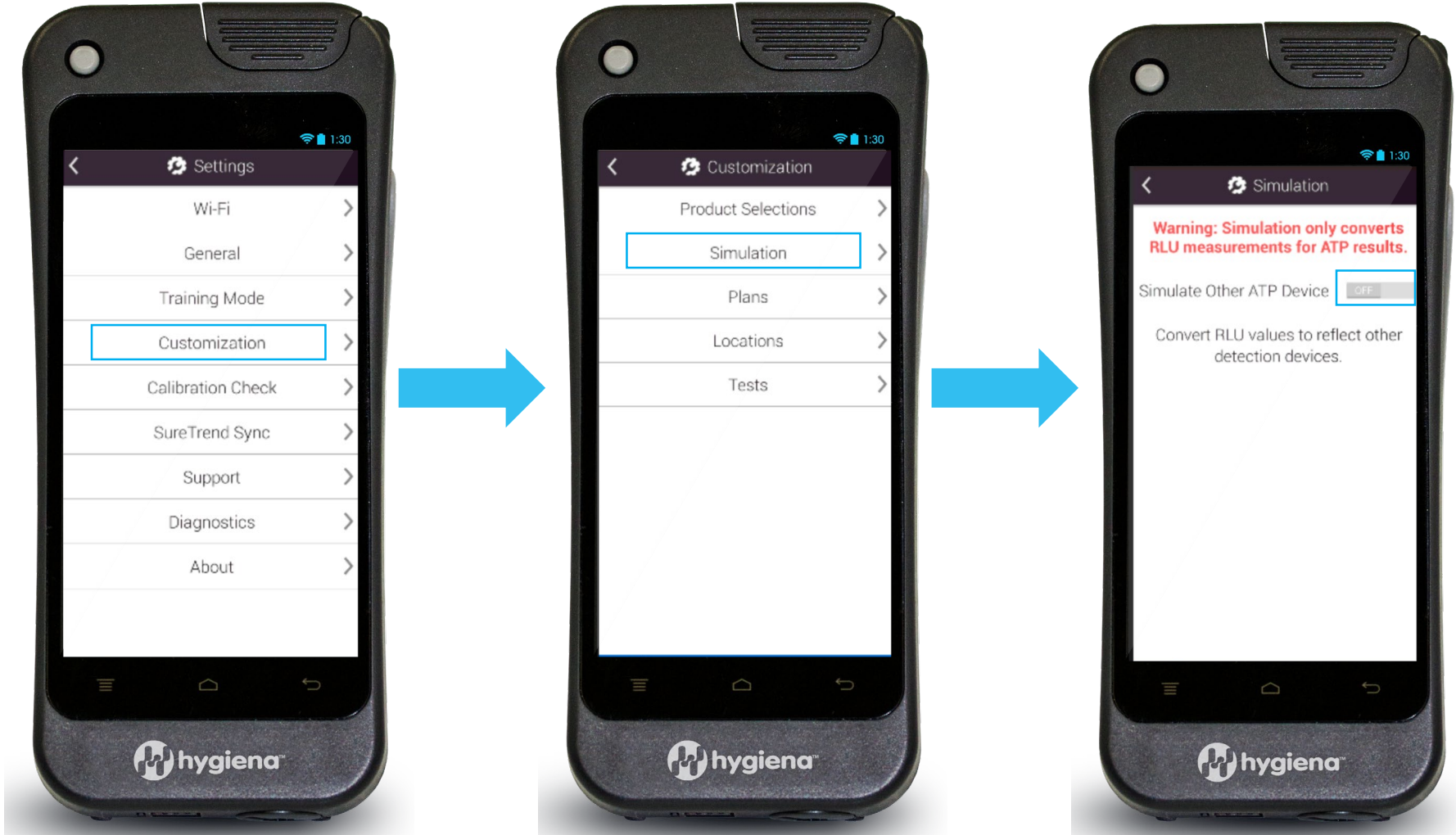




RLU Conversion

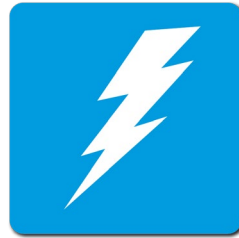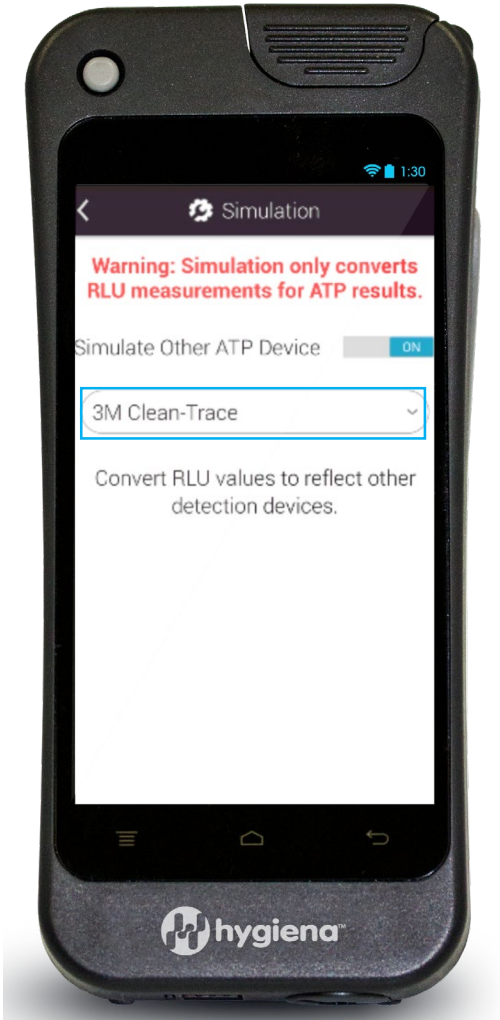

Run Test

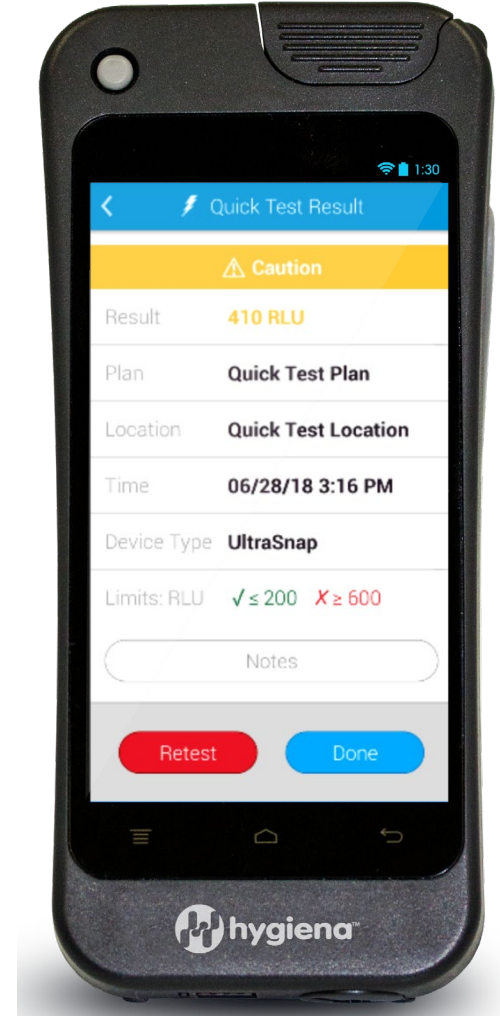

EnSURE™ Touch



Remote Support




- Built-in remote support on the instrument
- Easy screen sharing with a support representative
 - Screen sharing can be used for training employees





Remote Support

EnSURE™ Touch

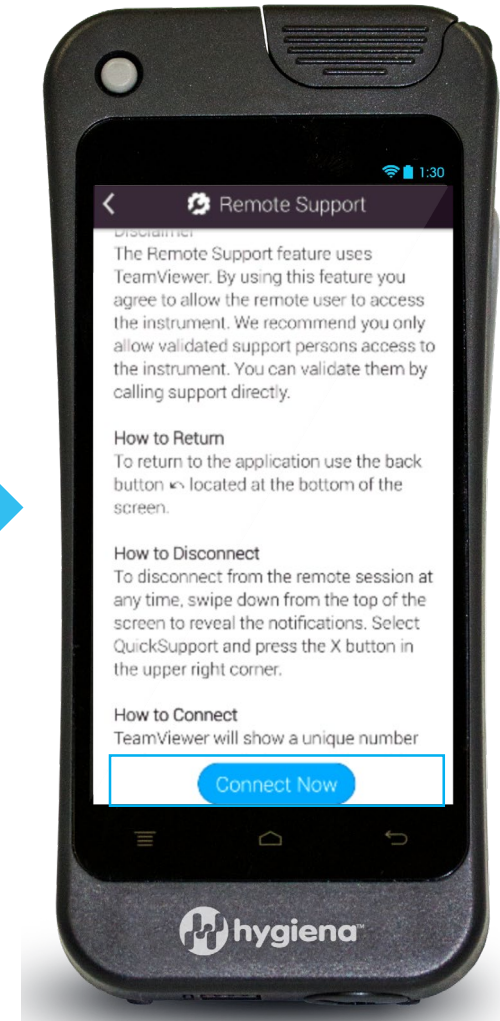
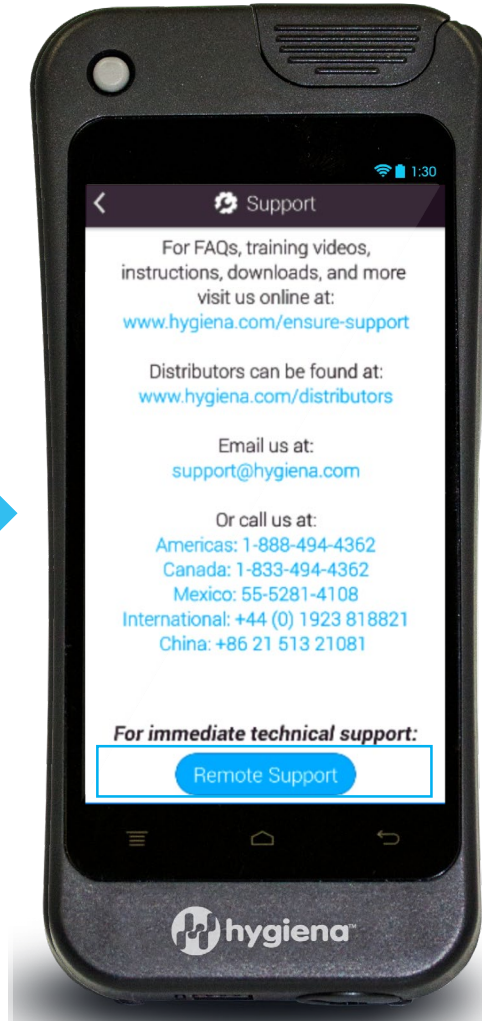
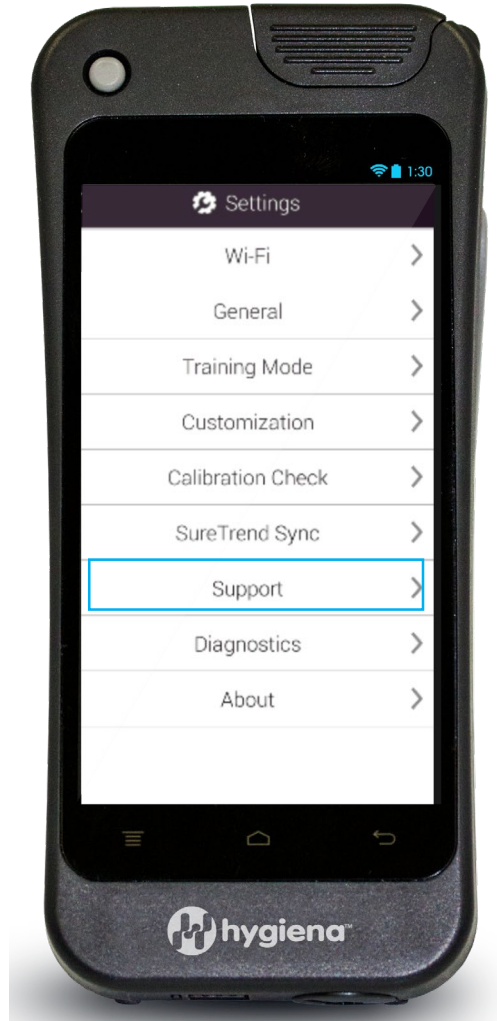





Remote Support Screen Share

EnSURE™ Touch

Customer's View

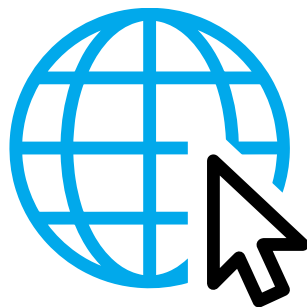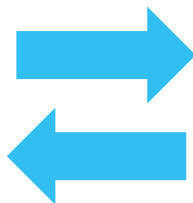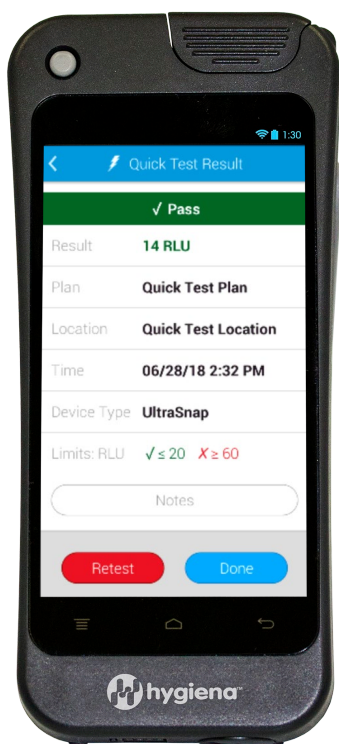


Support Representative's View

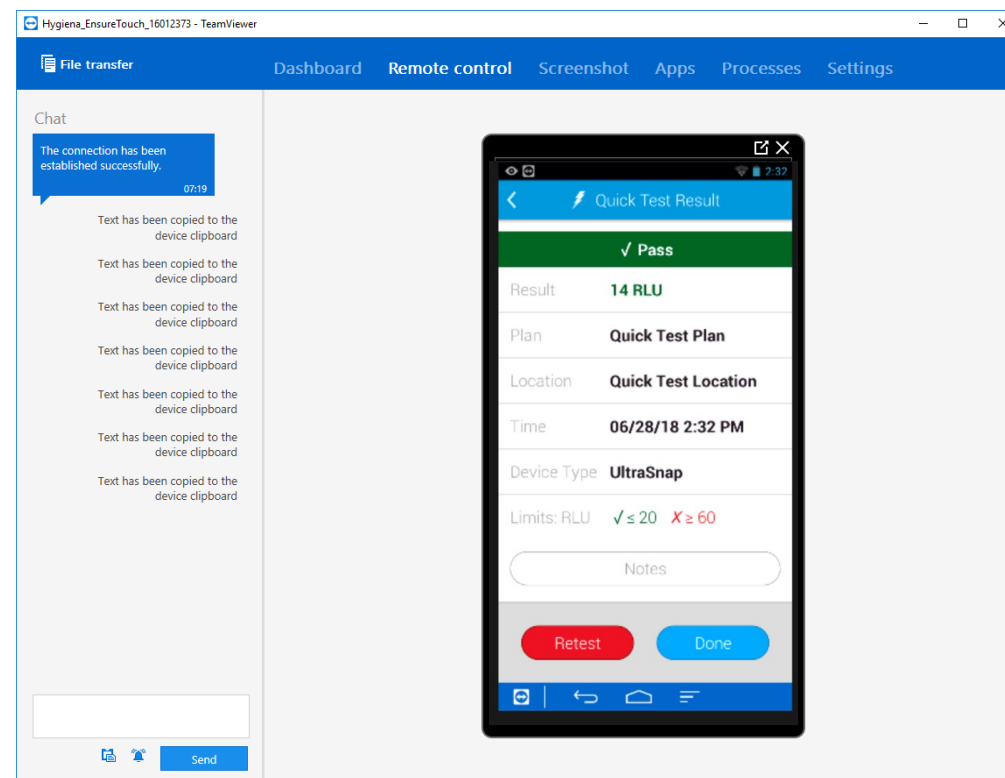



Welcome to SureTrend Cloud

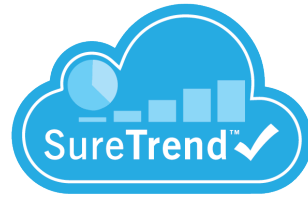

SureTrend[™] ✓ by  Hygiena[™]



Logout



Settings



Help

 Dashboard

 Results

 Reports

 Instrument

Welcome

SureTrend Cloud

Register
Instrument



Import Plans
and Locations



Migrate from
SureTrend[™]
V.4

Quick Support

Introduction to
EnSURE[™]
Touch

Video Tutorials



[Go To Dashboard](#)

☐ Don't show again

Bring your data to life with colorful & easy-to-use cloud software



SureTrend Cloud Dashboard



SureTrend™ by Hygiena™

[Logout](#) [Settings](#) [Help](#)


[Dashboard](#) [Results](#) [Reports](#) [Instrument](#)

Site: [Headquarters](#) Dashboard: [ATP - Fails By Plan](#) Period: [Custom](#) From: [1/1/2016, 12:00 AM](#) To: [6/4/2020, 11:59 PM](#)


ATP - Fails By Plan

Plan	Number Of Fails
Mill (All Fruit)	1143
	581
Quantarev	323
Mivap	304
Buhler Dryer	154

Pass Caution Fail By Plan: Mill (All Fruit)

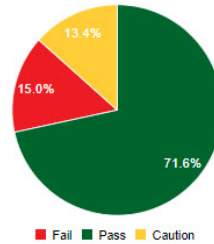

Pass Caution Fail Period

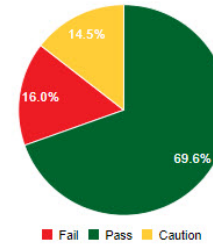


Fails For Location: Mill (All Fruit)

Location	Result Date	Rlu	Upper
Dumping Station Hop	12/14/2017 08:58:00 PM	377	30
Dumping Station Hop	12/14/2017 08:59:00 PM	220	30
Dumping Station Hop	12/14/2017 08:59:00 PM	222	30
Dumping Station Hop	12/15/2017 12:16:00 AM	62	30
Dumping Station Hop	01/07/2018 11:38:00 PM	103	30

1 2 3 4 5 ... 58

Pass Caution Fail By Plan: Mill (All Fruit)

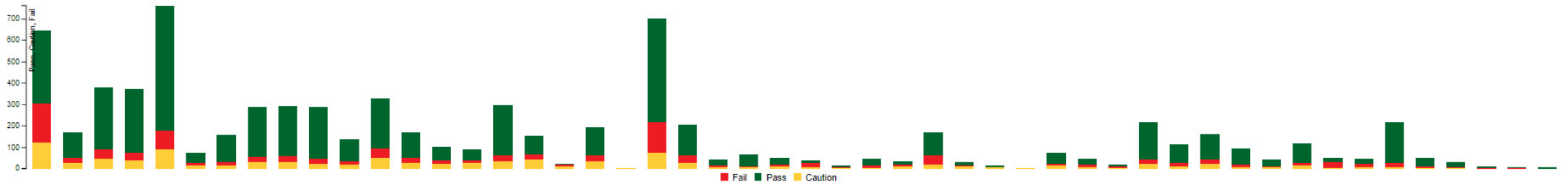


©2020 Hygiena, LLC

A bird's eye-view of sanitation practices across your organization



SureTrend Cloud Results

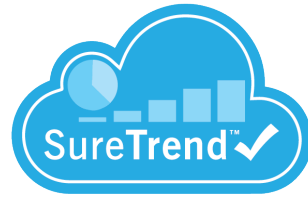

SureTrend™ by Hygiena

Logout Settings Help

Dashboard Results Reports Instrument

Type ATP

Period Custom

From 1/1/2016, 12:00 AM

To 6/1/2020, 11:59 PM

Drag a column header here to group by that column




	Result State	Result Date	Site Name	Location Name	Test Device	User Name		Rlu	Is Retest	Unit Name	Lower	Upper	Notes	Swab Temperature	A
✓	Pass	05/28/2020 12:26:...	Headquarters	Quick Test	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/21/2020 12:24:...	Headquarters	Computer keyboar...	UltraSnap			4	<input type="checkbox"/>	SEERGEI	100	100		18.3	
✓	Pass	05/13/2020 11:50:0...	Headquarters	Quick Test	UltraSnap	Renee		2	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/13/2020 11:49:0...	Headquarters	Quick Test	UltraSnap	Renee		2	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/13/2020 11:46:3...	Headquarters	Quick Test	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/12/2020 01:21:...	Headquarters	All Fill Chute	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	47.8125	60		0	
✓	Pass	05/12/2020 01:17:...	Headquarters	1	UltraSnap	Renee		2	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/12/2020 01:12:...	Headquarters	Box Filler Belt	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	48	60		0	
⚠	Caution	05/12/2020 12:37:...	Headquarters	Belt 1	UltraSnap	Renee		283	<input type="checkbox"/>	Product Managem...	280	350		0	
⚠	Caution	05/12/2020 12:37:...	Headquarters	Belt 1	UltraSnap	Renee		283	<input type="checkbox"/>	Product Managem...	280	350		0	
⚠	Caution	05/12/2020 12:36:...	Headquarters	Belt 1	UltraSnap	Renee		285	<input type="checkbox"/>	Product Managem...	280	350		0	
✓	Pass	05/12/2020 12:33:...	Headquarters	Quick Test	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/12/2020 12:30:...	Headquarters	Quick Test	UltraSnap	Renee		0	<input type="checkbox"/>	Product Managem...	20	60		0	
✓	Pass	05/12/2020 12:09:...	Headquarters	Quick Test	UltraSnap	Renee		4	<input type="checkbox"/>	Product Managem...	20	60		0	
✗	Fail	05/12/2020 12:09:...	Headquarters	Quick Test	UltraSnap	Renee		282	<input type="checkbox"/>	Product Managem...	20	60		0	
✗	Fail	05/12/2020 12:08:...	Headquarters	Quick Test	UltraSnap	Renee		283	<input type="checkbox"/>	Product Managem...	20	60		0	

8 12 20

Page 1 of 1580 (18958 items) 1 2 3 4 5 ... 1580


Database of all test results from all EnSURE Touch in your organization

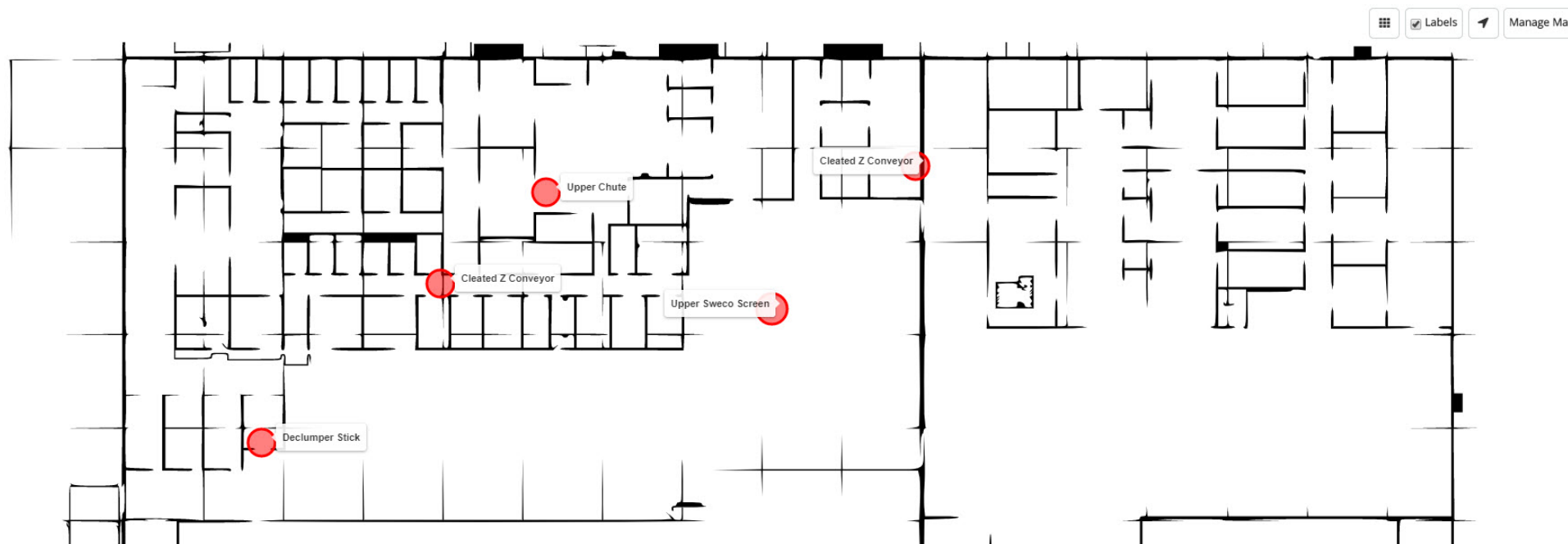

Track the origin of contamination with a visual indication of failed tests are on your floor plan



SureTrend Cloud Reports

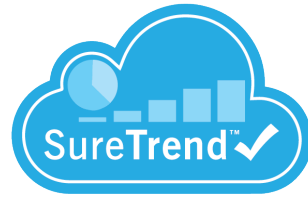

SureTrend™ by Hygiena

Logout Settings Help

Dashboard Results Reports Instrument

Reports

- Search
- Favorites
- Pass Caution Fail Summary
- Detail List
- Fails Summary
- Fails Not Retested
- Fails Retested
- Tests Performed
- Customized

Filters Options

Period: Custom From: 4/4/2016, 4:00 PM To: 4/11/2020, 3:59 PM

Plan: Select a value... Location: Select a value...


User: Select a value... Instrument: Select a value...

Result type: Pass ☐ Caution ☐ Fail ☐ Site: Select a value...

Retest: All results Deleted All results Apply Filter Reset Filter



PCF Summary by Day of Week

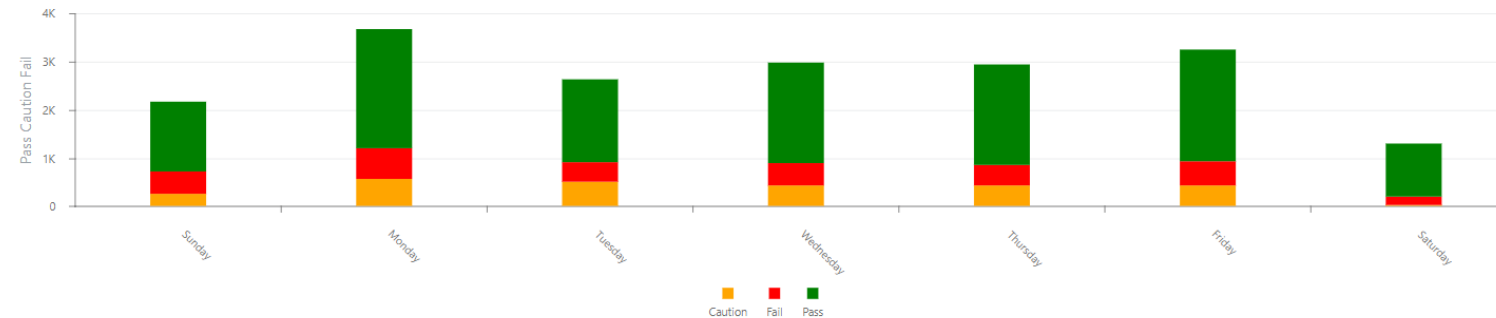


	Sunday	Monday	Tuesday	Wednesday	Thursday	Friday	Saturday
Caution	264	574	520	441	447	450	45
Fail	457	631	395	456	420	499	168
Pass	1443	2455	1724	2078	2067	2295	1102
Grand Total	2164	3660	2639	2975	2934	3244	1315

Pre-built & customized reports illustrate the success of your sanitation program



SureTrend Cloud Instruments

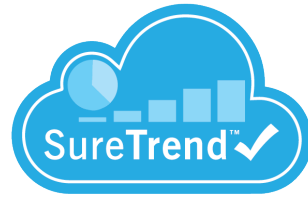
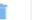

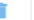
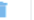







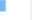






SureTrend[™] ✓ by 

 Logout
 Settings
 Help

 Dashboard  Results  Reports  Instrument

Instruments

	Serial No	Instrument Name	Site	Calibration Check	Model	Software	Last Sync
 	31325	Dev Team 1	Headquarters	Calibrated	EnsureTouch	1.39.12	22-Jul-2019 7:24 AM
 	33571	Product Management 2	Headquarters	Calibrated	EnsureTouch	1.55.7	24-Feb-2020 12:24 PM
 	30008	Product Management DE	Willmington, DE		EnsureTouch	1.48.3	10-Oct-2019 2:39 PM
 	31551	Product Management DE 2	Willmington, DE	Calibrated	EnsureTouch	1.41.10	22-Jul-2019 7:35 AM
 	32251	QA Lab 1	Seville, Spain		EnsureTouch	1.49.16	18-Oct-2019 2:36 PM
 	30843	QA Lab 2	Seville, Spain		EnsureTouch	1.42.9	25-Jun-2019 11:14 AM
 	31591	QC Lab	Mississauga, Canada	Calibrated	EnsureTouch	1.41.10	21-Jul-2019 2:53 PM
 	31040	R&D Lab 1	Watford, UK		EnsureTouch	1.48.3	18-Oct-2019 3:43 PM
 	30246	R&D Lab 2	Watford, UK	Calibrated	EnsureTouch	1.43.6	22-Jul-2019 7:23 AM
 	31690	Sales Team 1	Mexico City, Mexico		EnsureTouch	1.42.10	18-Oct-2019 1:17 PM
 	33536	Sales Team 2	Mexico City, Mexico	Calibrated	EnsureTouch	1.55.7	24-Feb-2020 12:32 PM

©2020 Hygiena, LLC

Manage & program your EnSURE Touch from anywhere in the world