EnSURE[®] Touch

Product Safety
Information

General

For technical assistance, please contact the Hygiena® location closest to your current region. For more information and technical support visit <https://help.hygiena.com>

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International:

Hygiena International
8, Woodshots Meadow
Watford, Hertfordshire
WD18 8YU, UK
Phone: +44 (0)1923 818821
Fax: +44 (0)1923 818825

Hygiena (Shanghai) Trading Co., Ltd.
Rm.7K, No.518, Shangcheng Rd.
Pudong New District
Shanghai, China 200120
Mobile: +86 13918834317
Phone: +86 21 51321081, 51321077/8/9
Fax: +86 21 51321080

Hygiena Diagnóstica España S.L.
P.I. Parque Plata – Calle Calzada
Romana,
40 · 41900 - Camas (Sevilla)
Phone: +34 955 983 215

Equipment Operation

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired. Use only as directed in the instructions located at <https://help.hygiena.com>

Equipment Maintenance and Service

All EnSURE® Touch components are supplied only by Hygiena or Hygiena's authorized agents and distributors (including, but not limited to, battery, power adapters, calibration devices, swabs). Only use Hygiena approved and distributed Lithium-Ion replacement batteries [MIS0096]. Use of any other battery may cause damage to the device. Please refer to your local Hygiena Sales Representative or <https://help.hygiena.com> for help placing your replacement order.

Reboot Necessity & Instructions for Exposure to High Frequency Radiation and Electrostatic Discharge

This device is susceptible to Electrostatic Discharge (ESD). In the presence of ESD an operator may experience display degradation and/or loss of touch screen function. Exposure to high frequency radiation may cause the EnSURE Touch instrument to perform unexpectedly and outside of normal operating function. If this occurs, either remove the source of radiation, or remove the instrument from the source of radiation. Then perform a hard reboot of the instrument by holding the power button at top left for three (3) seconds until the instrument restarts. If the instrument continues to exhibit unexpected, or outside of normal operating function, please notify your local Hygiena representative.

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Hygiena Mexico, S.A. de C.V.